



SCILOGEX D160/D500 Homogenizers





Tel: (860) 828-5614 Fax: (860) 828-5389 info@scilogex.com www.SCILOGEX.com





The Scilogex D160 Homogenizer can rapidly homogenize, disperse, and emulsify samples in 0.1 - 50 ml of liquid or 1-250ml depending on the generating probe. During operation, the suspended material is drawn into the core of the homogenizer by a rotor turning at up to 30,000 rpm. The material is repeatedly cycled through narrow slits in the probe where it is rapidly sheared and disintegrated by high shear. Complete homogenization of tissues (muscle, liver, breast tissue, etc.) is usually achieved in a few seconds. Little, if any, heat is produced during the process.

Features:

Use for mixing, emulsification, shredding, or chopping.
Practical answer for small samples.
Easy to hold and light weight for one hand operation.
Generators are made of 316 L Stainless Steel with Teflon bushings.
Autoclavable stainless steel shafts. Inert to aggressive solutions.
All shafts are interchangeable shafts and are easily removed for cleanup.
High speed motor. Speed control from 8,000 to 30,000 rpm.
Suitable for solvents & abrasive substances.

Specifications:

Working Volume (ml): 0.1 - 250ml

Operating Environment: 0-40°C, 95% rel. humidity
Motor type: Universal motor, 130 Watt

Voltage / Frequency: 110V/60Hz

Certified standards: 89/336EEC,72/023/EEC, EMC acc. IEC/EN 61000-6-2 / 3,

Safety acc. IEC/EN 61010-2-51

Speed range (rpm): 8,000 - 30,000 Warranty: 2 years

Speed control: 6 position control dial Noise emission: < 54 dB(A) at 30,000 rpm

pH Range: 2-13

Dimensions (mm): 250L x 250W x 400H

Weight (kg) 1.6

Fineness, (µm) Suspensions & Emulsion 10-15 & 1-10

Ordering Information:

Cat No.

Description

SCILOGEX D160 Homogenizer Package, DS-160/5 5mm dia. generator, flat head-open slot, 50mm immerse length, H-370 stand, for solid/liquid media, volumes 0.1-50ml, 110V, 60Hz

SCILOGEX D160 Homogenizer Package, DS-160/10 10mm dia. generator, flat head-open slot, 115mm immerse length, H-370 stand, for solid/liquid media, volumes 1- 250ml, 110V, 60Hz

85010301 SCILOGEX D160 Homogenizer Drive Motor, 110V/60Hz

85010401 SCILOGEX H-370 Homogenizer Stand

Accessories:

18900113 DS-160/5 5mm dia. generator, flat head-open slot, 50mm immerse length for D-160,

for solid/liquid media, volumes 0.1-50ml

18900114 DS-160/10 10mm dia. generator, flat head-open slot, 115mm immerse length for D-160,

for solid/liquid media, volumes 1- 250ml









The Scilogex D500 Homogenizer has a maximum speed of 29,000 rpm rate, and educes processing time significantly. Smooth motor drive makes working, very safe and comfortable. An overload protection feature increases the life span of the motor. Smooth acceleration also prevents unnecessary spills. Flexible configuration of different generators and rotors make the D500 highly adaptive to most applications. The generator shaft is constructed of quality 316L stainless steel. Reliable results are assured with extremely high precision and speed. Assembly and disassembly of rotors requires no special tools.

Specifications:

Sample Volume Range: 1 to 50ml or 10 to 5000ml Universal Motor: 110V/60 Hz, 500 watts

Variable Speed: 10,000 to 29,000 rpm, incremented by scale

Weight: 1.6 Kg (drive only)
Operating Environment: 0-40°C, 95% rel. humidity
Motor type: Universal motor, 500 Watt

Voltage / Frequency: 110V/60Hz

Certified standards: 89/336EEC,72/023/EEC, EMC acc. IEC/EN 61000-6-2 / 3,

Safety acc. IEC/EN 61010-2-51

Speed range (rpm): 10,000 - 29,000

Warranty: 2 years

Speed control: 6 position control dial Noise emission: < 72 dB(A) at 30,000 rpm

pH Range: 2-13

Dimensions (mm): 250L x 250W x 700H Weight (kg) 9.0 Kg (with stand)

Fineness, (µm) Suspensions & Emulsion 10-15 & 1-10

Ordering Information:

Cat No.	Description

85020101 SCILOGEX D500 Homogenizer Package, DS-500/1 20mm dia. generator, with S20C/SR20 flat head-open

slot coarse generator, 215mm immerse length, for solid/liquid media, volumes 10-5000ml and with

H-600 stand, 110V, 60Hz

85020201 SCILOGEX D500 Homogenizer Package, DS-500/2 20mm dia. generator, with S20F/ER20 flat head-open

slot fine generator, 215mm immerse length, for water in oil or oil in water emulsions, volumes

10-50000ml and with H-600 stand 110V, 60Hz

85020301 SCILOGEX D500 Homogenizer Drive Motor, 110V/60Hz

85020401 SCILOGEX H-600 Homogenizer Stand

Accessories:

18900115 DS-500/1 20mm dia. generator, with S20C/SR20 flat head-open slot coarse generator, 215 immerse length

for solid/liquid media, volumes 10-5000ml

18900116 DS-500/2 20mm dia. generator, with S20F/ER20 flat head-open slot fine generator, 215 immerse length

for Water in Oil or Oil in Water emulsions, volumes 10-5000ml

18900117 DS-500/5 generator 5mm dia. generator, flat head-open slot, 60mm immerse length for solid/liquid

media, volumes 1-50ml

18900118 S20C 20mm open slotted coarse generator 18900119 S20F 20mm open slotted fine generator 18900120 S30C 30mm open slotted coarse generator 18900121 S30F 30mm open slotted fine generator 18900122 S40C 40mm open slotted coarse generator

18900123 SR20 20mm standard rotor SR30 30mm standard rotor 18900124 CR20 20mm knife rotor 18900125 CR30 30mm knife rotor 18900126 18900127 MR20 20mm mixing rotor 18900128 MR30 30mm mixing rotor 18900129 ER20 20mm emulsifying rotor 18900130 ER30 20mm emulsifying rotor





18900113









Generators

18900115











Rotors











18900125



18900126



18900127

Rotors







18900129



18900130

